

# Chromium(VI)

## Test kit for performing colorimetric tests on chromium(VI) in surface water and sewage

### Method:

In an acidic medium chromate ions react with diphenylcarbazide to form a red-violet dye. First chromium(VI) oxidizes diphenylcarbazide to diphenylcarbazone, being itself reduced to chromium(III). Combined with the enol form of the carbazone, these chromium(III) ions form the intensively coloured complex.

### Measurement range:

0.02 - 0.50 mg/l Cr(VI)

### Contents of test kit (\*refill pack):

sufficient for 140 tests

30 ml Cr-1\*

25 ml Cr-2\*

2 screw-plug measuring glasses

1 slide comparator

1 colour chart

1 plastic syringe 5 ml

1 instructions for use\*

### Hazard warning:

Cr-1 contains *ortho*-phosphoric acid 68%. **Causes burns.** In case of contact with eyes, rinse immediately with plenty of water and seek medical advice. For further information, please ask for a safety data sheet.

### Instructions for use:

also refer to the pictogram on the back of the colour chart

1. Pour a 5 ml water sample into each of the measuring glasses using the plastic syringe.  
Place a measuring glass on position A in the comparator.

### Only add the reagent to measuring glass B.

2. Add **5 drops of Cr-1**. Seal the glass and mix.
3. Add **5 drops of Cr-2**. Seal the glass and mix.
4. Open the glass after **3 min** and place it on position B in the comparator.
5. Slide the comparator until the colours match in the inspection hole on top. Check the measurement reading in the recess on the comparator reed. Mid-values can be estimated.
6. After use, rinse out both measuring glasses thoroughly and seal them.

The method can also be applied for the analysis of sea water.

The reagents can also be used for the **photometric evaluation** with photometer PF-11.

### Disposing of the samples:

The used analysis specimens can be flushed down the drain with tap water and channelled off to the local sewage treatment works.

### Interferences:

Only chromium(VI) can be determined with this test kit. Chromium(III) must first be oxidized (see below determination of *total* chromium). Larger quantities of heavy metal ions interfere.

### Conversion table:

mg/l Cr(VI)	mg/l CrO <sub>4</sub> <sup>2-</sup>
0.02	0.04
0.05	0.11
0.10	0.22
0.15	0.33
0.20	0.45
0.30	0.67
0.40	0.89
0.50	1.12

### Determination of *total* chromium:

Add to 20 ml test sample 1 ml sulphuric acid 96% and 0.5 g potassium peroxodisulphate and boil for 2 h. After cooling add approximately 10 ml of distilled water and carefully adjust to pH 1-3 with 5 ml of sodium hydroxide solution 20%. Then fill up to 50 ml and determine the chromium concentration as described above. Multiply the read-off value with **2.5**.

### Note:

For the determination of water-soluble chromium(VI) in cement contact MACHEREY-NAGEL for special working instructions.

### Storage:

Store the test kit in a cool (< 25 °C) and dry place.